**ritter**  
MEDICAL

# Liquid Handling

Ripette® pro,  
Ritips® evolution  
and more.

### Liquid-Handling-Systeme

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### Liquid handling systems

In order to make high precision dosing easier and faster, Ritter Medical offers repetitive dispensers for manual use which allow a repeatedly dispensing of a specific aliquot. For the dosing of different aliquots, Ritter offers a wide range of precisely engineered consumables which are available also in a bioclean® (sterile) grade. Our target is to offer you one dosing system which guarantees reliable and precise results for many applications.

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## From a first product idea to just-in-time delivery

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With an overall target of maintaining healthy life, Ritter Medical supports research as well as diagnostics with new product innovations. From the first product idea to the development on modern CAD systems and in-house tool and mould production, Ritter creates new products in order to offer the best solution for your applications.

All project processes are combined under one umbrella, i.e. injection moulding, quality assurance as well as logistics and distribution. In more than 70 countries, millions of products of Ritter Medical are used in clinical and biotechnological applications on a daily basis. Furthermore, our products are used in numerous applications within the area of personalized medicine as well as for applications which help to make routine processes within a laboratory safer and easier. It is our target to always offer you an optimal value for money at the same time!

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Ripette® pro is ergonomically the best choice by offering a high flexibility and handy features. It is furthermore designed and constructed for mechanical dispensing of highly precise dosage volumes systematically. Due to its mechanical construction, the Ripette® pro becomes low maintenance and requires no additional resources such as a power supply for charging batteries.

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The Ripette® pro is optimized for dispensing of contaminated fluids and volatile liquids. Using Ritips® evolution and other compatible, the Ripette® pro is reliably protected against contamination by aerosols ascending from the tip. Also on a long-term it can be securely used even with highly fluid samples.

The Ripette® pro offers a very wide choice of dosing volumes, allowing quick tip changes between 12 tip sizes. By using the 10-stages adjustment ring, up to 120 different program steps can be chosen in order to dispense volumes from 1 µl up to 5,500 µl.

The ergonomic shape ensures a user-friendly and easy operation. The low weight (87g) combined with a dispensing button located on top of the device allow a fast and easy dispensing even of long series.

## Advantages

- **Great precision:** The Ripette® pro's vision panel clearly shows the currently used setting as well as the dosing volume for the particular tip size.
- **High compatibility:** Ripette® pro is precisely aligned with Ritips® and Ritips® evolution from Ritter but can also be used with Brand's PD-tips and Eppendorf's Combitips advanced®.
- **Robust:** By using only high-quality materials for the production of Ripette® pro, the dispenser offers an excellent chemical resistance and a long service life.
- **Tested quality:** Every device is individually tested for its functionality and precision after production. Only if the requirements of DIN EN ISO 8655 have been fulfilled, the Ripette® pro will get an inspection certificate with its serial number.

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The Ripette® has been designed for dispensing precise volumes in long series. At the same time a high value has been set to make it ultra-robust and mechanically reliable by using only a minimum of moving parts. Due to its light weight (105g) it is well-suitable for long dispensing series. The ergonomic design allows a single-handed volume selection, loading and dispensing. The user can quickly select the ideal setting for a particular application, choosing either the highest precision or the maximum number of repetitions of the desired dosage volume.

When used in combination with Ritips® and Ritips® evolution, the attached dosage guide allows the user to define the best tip for his application at a glance.

Each device is tested for its functionality and precision according to DIN EN ISO 8655 before a serial number is applied which ensures a twelve-month warranty.

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Ritips® are available in eight classic sizes which include also the 'in-between' sizes 0.05 ml, 1.25 ml and 12.5 ml. By using the direct displacement method, Ritips® work well for infectious and volatile liquids, as the fluid never will get in touch with the instrument which avoids any contamination.

For sensitive applications or analysis of body fluids, Ritips® are available also as a bioclean® grade (single sealed and sterilized). Automated individual tests of each single tip guarantee the tightness with water, which is documented in the lot-specific records.

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Ritips® evolution are engineered to provide universal syringes in line with the Ripette® pro as well as with most other dosing systems. Ritips® evolution are compatible with several steppers for different applications. It is our target to offer you just one dosing system which works reliably and precisely for all your applications.

The Ritips® evolution consist of nine tip sizes for repeated dispensing volumes, ranging from 1µl to 50ml. Every single tip is subject to automated, individual testing to ensure 100% reliable and repeatable results.

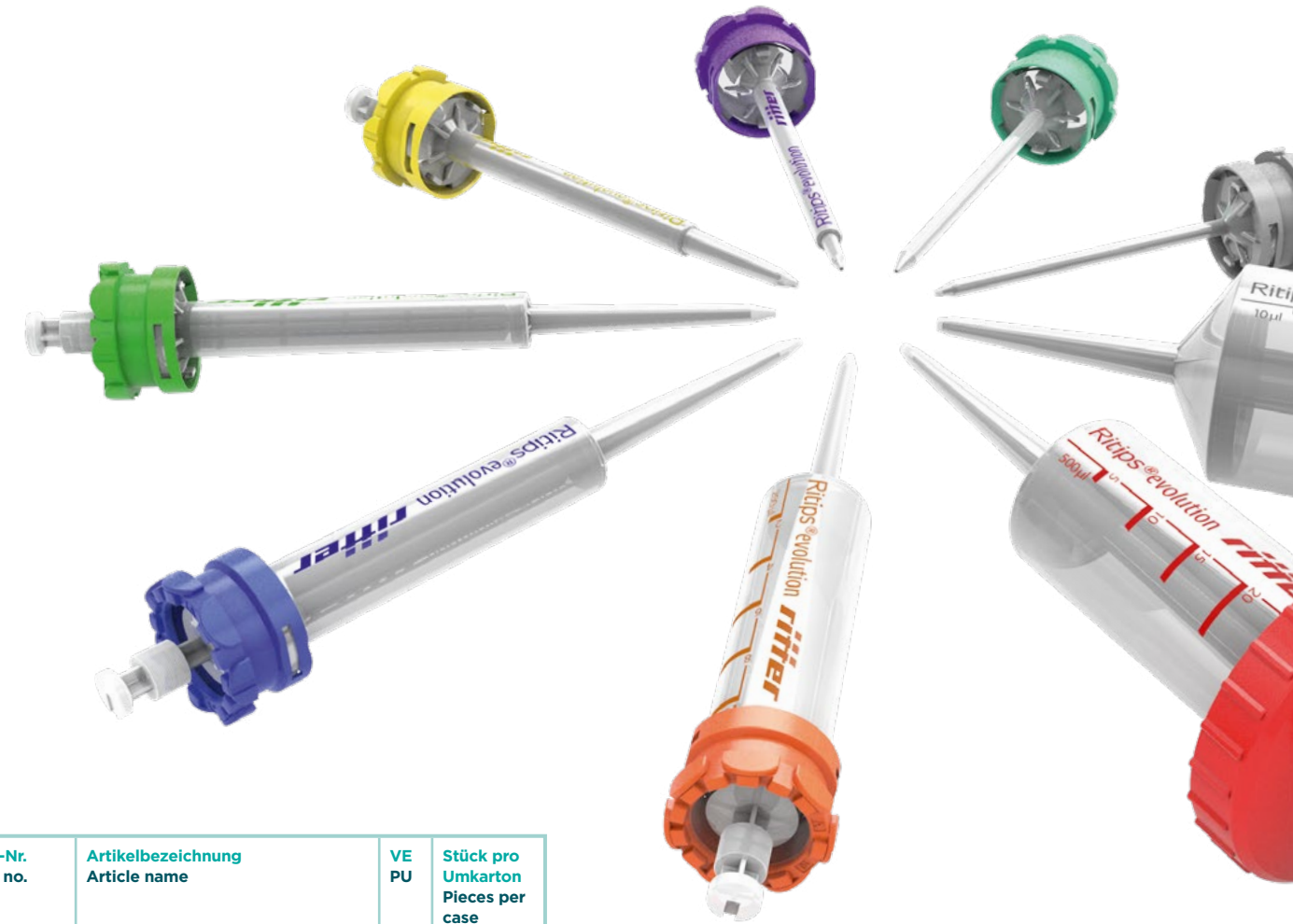
Ritips® evolution are available in standard quality in packs of 25 and 100 pieces, respectively, as well as in bioclean® quality: sterile and single-wrapped. This can be confirmed by independent laboratories which certify that the tips are sterile and free from endotoxins, ATP, RNase and DNA impurities. Certificates for each batch are available on request, with lot number recorded on each box for full traceability.

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The mini bottle dispenser is used for quick and easy dispensing of liquids in volumes of either 500 µl or 1000 µl, its standard thread size is 1 inch (26 mm). The output volume can be selected by twisting an adjustment wheel and then operating the centrally mounted plunger.

After refilling, the bottle dispenser can be vented through repeated pumping of the piston. The bottle is then ready for repeated use.

The bottle dispenser has no metal valves and is made from shatterproof polypropylene with very high chemical resistance. Cleaning can be done by a damp cloth, using soapy water solution if necessary.

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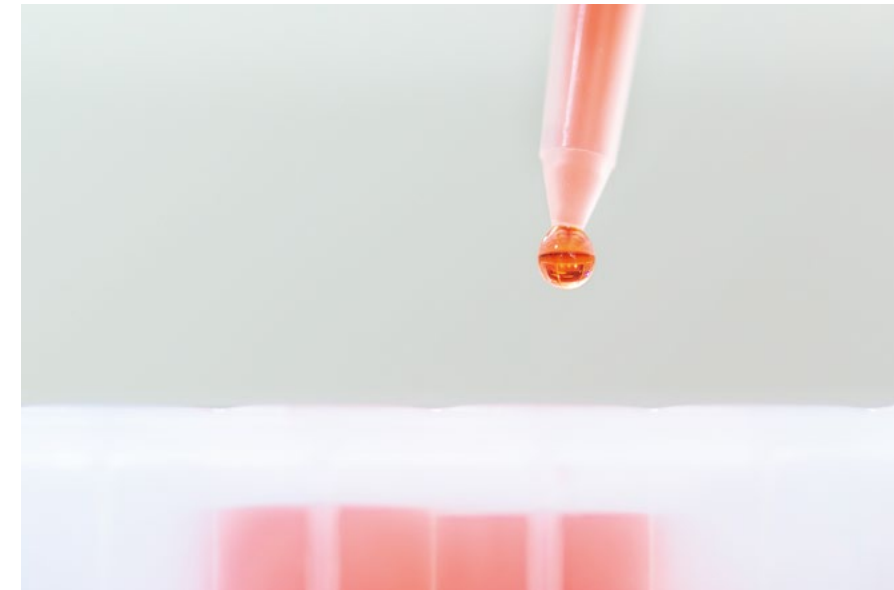
According to DIN EN ISO 8655, dispensers have to be controlled minimum once per year. Therefore, you can send your devices including a calibration order to Ritter Medical. We are glad to take care about the inspection of your or your customer's dispenser. Nevertheless, we will leave the decision about the inspection of your devices up to you. No matter if you would like to make the calibration of your dispenser in-house or by yourself, please follow the detailed instruction of DIN EN ISO 8655-1/7.3. The enclosed chart provides the necessary technical data for a successful calibration.

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The data for Ritips® 0.05 ml in the chart are based on a measurement with a 10 µl pipette tip attached. Measurement according to DIN EN ISO 8655 with Ripette® pro. Liquid: bidistilled water. Reference temperature: approx. 20 °C. The Ripettes are suited for a temperature range from +15 °C to +40 °C.

	Dispenservolumen in µl Dispensed volume in µl			Systematische Messabweichung in % Measurement deviation determined by the system in %			Zufällige Messabweichung in % Incidental measurement deviation in %		
	2% V <sub>nom</sub>	6% V <sub>nom</sub>	10% V <sub>nom</sub>	2% V <sub>nom</sub>	6% V <sub>nom</sub>	10% V <sub>nom</sub>	2% V <sub>nom</sub>	6% V <sub>nom</sub>	10% V <sub>nom</sub>
Tipgröße Tip size									
0.05 ml	1	3	5	± 6	± 4	± 4	≤ 9.0	≤ 5.0	≤ 4.0
0.10 ml	2	6	10	± 2.5	± 2	± 1.5	≤ 4.5	≤ 4.0	≤ 2.0
0.20 ml	4	12	20	± 1.6	± 1.6	± 1	≤ 1.9	≤ 1.2	≤ 1.0
0.50 ml	10	30	50	± 1	± 1	± 1	≤ 1.5	≤ 1.2	≤ 0.8
1.00 ml	20	60	100	± 1	± 1	± 1	≤ 1.3	≤ 1.2	≤ 0.9
1.25 ml	25	75	125	± 0.9	± 0.9	± 0.9	≤ 0.9	± 0.7	≤ 0.6
2.50 ml	50	150	250	± 0.7	± 0.7	± 0.7	≤ 0.9	≤ 0.7	≤ 0.6
5.00 ml	100	300	500	± 0.3	± 0.3	± 0.3	≤ 0.8	≤ 0.6	≤ 0.4
10.00 ml	200	600	1000	± 0.3	± 0.3	± 0.3	≤ 0.8	≤ 0.6	≤ 0.4
12.50 ml	250	750	1250	± 0.3	± 0.3	± 0.3	≤ 0.5	≤ 0.5	≤ 0.4
25.00 ml	500	1500	2500	± 0.5	± 0.5	± 0.5	≤ 0.6	≤ 0.4	≤ 0.3
50.00 ml	1000	3000	5000	± 0.5	± 0.5	± 0.5	≤ 0.6	≤ 0.5	≤ 0.3

# Qualitätsmanagement Quality Management

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In more than 70 countries, millions of plastic products of Ritter are used in clinical and biotechnological applications on a daily basis. In order to guarantee a high quality of products, Ritter has been certified according to ISO 9001. For the production of medical and IVD-products Ritter is also certified in line with the ISO 13485:2016.

For the production of IVD products, Ritter applies only certified and virgin raw materials. Sustainability and performance according to the EnMS ISO 50001:2011 are significant for a responsible manufacturing process.

## Certificate

Standard **ISO 9001:2015**  
Certificate Registr. No. **01 100 120473**

Certificate Holder:  
**ritter**  
Ritter GmbH  
Kaufbeurer Str. 55  
86830 Schwabmünchen  
Germany

Scope: Design, development, production and distribution of plastic components, particularly of cartridges, medical articles, laboratory products and building materials

Proof has been furnished by means of an audit that the requirements of ISO 9001:2015 are met.

Validity: The certificate is valid from 2019-01-17 until 2022-01-16. First certification 2013

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www.tuv.com

## Certificate

The Certification Body of  
TÜV Rheinland LGA Products GmbH

hereby certifies that the organization  
**Ritter GmbH**  
Kaufbeurer Str. 55  
86830 Schwabmünchen  
Deutschland

has established and applies a quality management system for medical devices for the following scope:  
**Design and development, manufacture and distribution of dialysis cartridges as well as accessories for In-Vitro Diagnostics (other IVD from plastic) and plastic injection moulding articles for medical devices (non-implants)**

Proof has been furnished that the requirements specified in  
**EN ISO 13485:2016**  
are fulfilled. The quality management system is subject to yearly surveillance.

Effective Date: 2019-03-22  
Certificate Registration No.: SX 80137032 0001  
An audit was performed. Report No.: 21229000 008  
This Certificate is valid until: 2019-11-29

Certification Body

Date 2019-03-22

**TÜV Rheinland LGA Products GmbH - Tillystraße 2 - 90431 Nürnberg**  
Tel.: +49 231 858-1271 Fax: +49 231 908-3929 e-mail: cert-verify@tuv.com http://www.tuv.com

## MANAGEMENT SYSTEM CERTIFICATE

Certificate No. 134459-2013-AE-GER-DAAS Initial certification date: 28. June 2013 Valid: 28. June 2016 - 27. June 2019

This is to certify that the management system of  
**Ritter GmbH**  
Kaufbeurer Str. 55, 86830 Schwabmünchen, Germany

has been found to conform to the Energy Management System standard:  
**ISO 50001:2011**

This certificate is valid for the following scope:  
**Design, development, production and distribution of plastic components, particularly of cartridges, medical articles, laboratory products and building materials**

Place and date:  
**Essen, 30. May 2018**

For the issuing office:  
**DNV GL - Business Assurance**  
 Schöteninghof 14, 45328 Essen, Germany  
  
**Thomas Beck**  
 Technical Manager

Lack of fulfillment of conditions as set out in the Certification Agreement may render this Certificate invalid.  
Tel.: +49 201 724-222 www.dnv.com

**Ritter GmbH**

Kaufbeurer Str. 55  
86830 Schwabmünchen  
Germany

phone: +49 8232 5003-45  
[www.ritter-medical.de](http://www.ritter-medical.de)

*ritter*

93002-0164 | V19